

OMERO search

Step-by-step guide: How can I **search** for data of specifics user?

1. In the top bar of the OMERO web page you will find a "Search" field on the right side. There you can enter a full name to search for.
2. You can filter the results for specific groups, object types or other restrictions.
3. Be aware that using wildcards (see yellow box) is possible and might be needed, in case you are not searching for an exact name.

The screenshot shows the OMERO webclient interface with a search query of "Anna Hamacher". The search results are displayed in a table with columns: Type, Name, Acquired, Imported, and Group. The results list various datasets and projects associated with the user Anna Hamacher.

Search Interface Details:


- Search Bar:** Located at the top right, containing the text "Anna Hamacher".
- Filter Results:** A dropdown menu showing "More than 520 results".
- Search Hints:** A yellow box on the left provides information about wildcards: "Single character wildcard: ?" and "Multiple character wildcard: *". It also states "Results will contain both terms. E.g: GFP AND H2B".
- Search Field:** A text input field containing "Anna Hamacher".
- Restrict by Field:** Checkboxes for Name, Description, and Annotations.
- Search for:** Checkboxes for Images, Datasets, Projects, Wells, Plates, and Screens.
- Scope:** A dropdown menu set to "All Groups".
- Data owned by:** A dropdown menu set to "Anna Hamacher".
- Date:** A dropdown menu set to "Import date".
- Search Button:** A blue button labeled "Search".

Search Results Table:


Type	Name	Acquired	Imported	Group
Image	CAI_ITIC_2021_Anna	2021-04-01 09:47:56		CAI_CAI
Image	2021-Feb_LSM880_Airyscan	2021-03-03 11:07:03		CAI_Membra
Image	Demo_Wiki	2021-05-10 12:26:43		CAI_CAI
Image	FM464-FX	2021-03-03 11:07:24		CAI_Membra
Image	Cellmask_Phototox_Blebbing	2021-03-03 11:07:37		CAI_Membra
Image	Cellmask_after4days_Phototox_Blebbing	2021-03-03 11:07:13		CAI_Membra
Image	Fixed_Tests	2021-04-01 09:49:34		CAI_CAI
Image	Live_Pretests_MbxA	2021-04-01 09:49:26		CAI_CAI
Image	Live_MbxA_old_3days	2021-03-24 14:12:11		CAI_CAI
Image	Live_Calcein	2021-04-01 10:00:28		CAI_CAI
Image	Live_CalciumSensor	2021-04-01 09:49:16		CAI_CAI



If you are actually not sure who the owner of the data is, you are interested in, you can browse the Data of a group by using the "All Members" navigation.

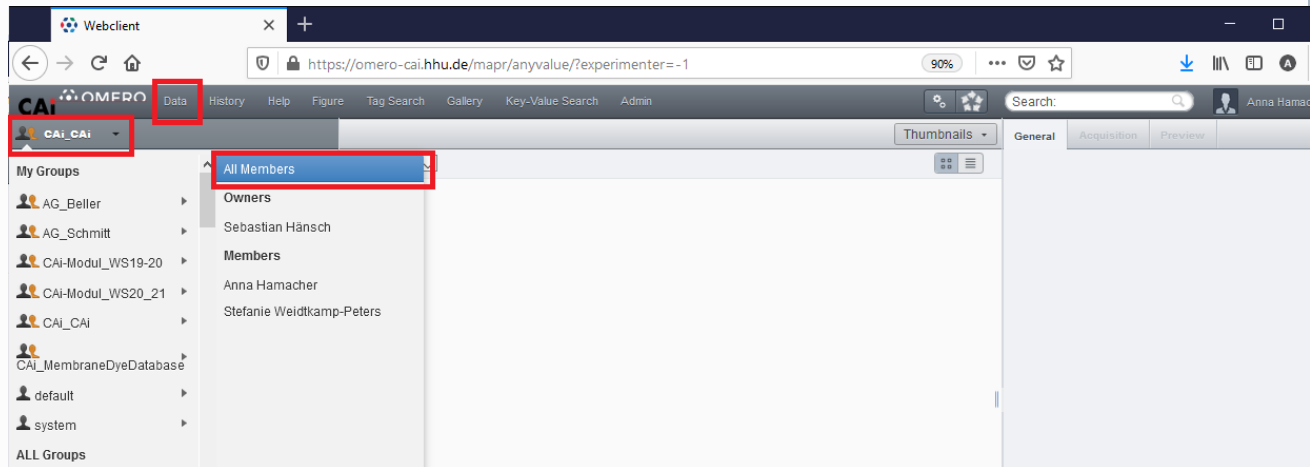

Alternatively you can use the tag search or key-value search if these annotations are maintained.

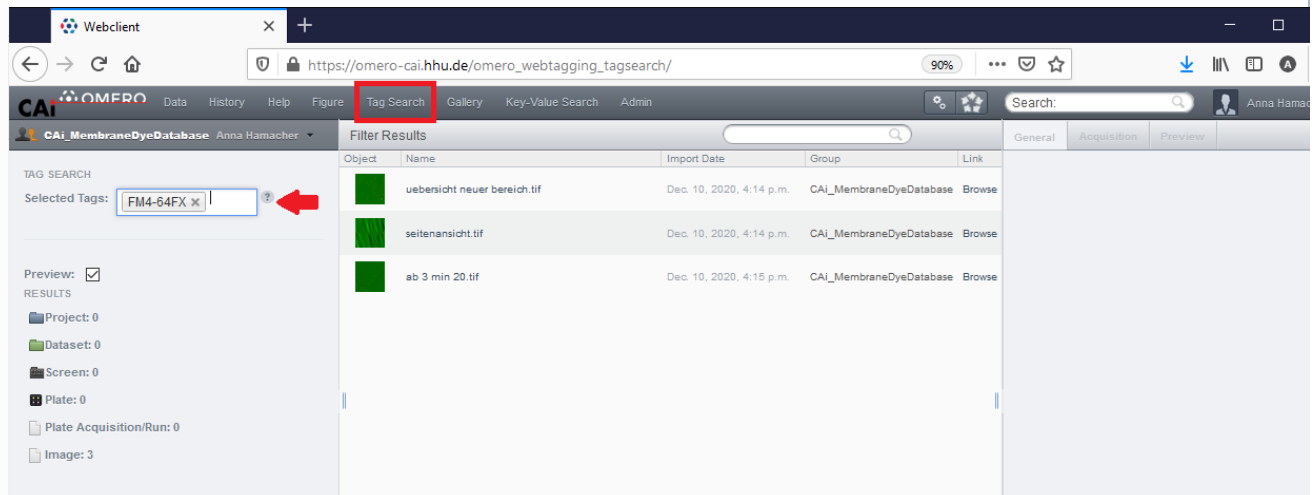

Related articles

- [OMERO download](#)
- [OMERO upload](#)
- [OMERO search](#)
- [OMERO web link](#)
- [OMERO data sharing](#)