OMERO group owner

As a group owner you have additional permissions within your OMERO group. One important permission is the management of the users within your group.

Step-by-step guide: I'm a group owner, how can I modify the users in my group?

• Left-click on your profile (top-right within the OMERO web page) > "User Settings".

	😧 Webclient	× +		- 0
$\left(\leftarrow \right)$	ightarrow C $harrow$	🖲 🔺 https://omero-cai. hhu.de /webclient/	90% 🗹 🏠	👱 III\ 🗊 🔕
CA	Data History	Help Figure Tag Search Gallery Key-Value Search Admin	Search:	🔍 🚶 Anna Hama
2 c	Ai_MembraneDyeDatabase Anna	amacher *	Thumbnails - General Acquisitio	User settings
Explor	Tags Shares			Send Feedback
				- Logout
- 🗶 Ar	na Hamacher 2021-Feb_LSM880_Airyscan 3			
	Orphaned Images 1			

· Choose the tab "My Groups" and click on the "Edit" button related with the group you would like to modify.

🔅 My account	× +				- 0
← → ♂ ☆	🛛 🔒 https://omero-cai.hhu.de/	/webadmin/myaccount/edit/		90% 🗹 🏠	⊻ III\ 🗊 🔕
CALIFIC Data History	Help Figure Tag Search Galle	ery Key-Value Search Admin		Search:	🔍 🚶 Anna Hama
	User Wy Groups 3	Permissions Read-Annotate (wra-) Read-Annotate (wra-)	Description - For demo and documentation purposes	Edit	

- Left-click into the "Members" field will open a list of persons you can add to your group. By writing a part of the name into this field, the list of persons will be filtered.
- If you found the person you would like to add, click on it to add it to the "Members" field.
- Afterwards use "Save" to make your change permanent.

😧 Edit group 🛛 🗙	+		- 0
$\leftarrow \rightarrow$ C $\textcircled{0}$	https://omero-cai.hhu.de/webadmin/group_owner/edit/1358/	F 90% ··· 🗸	👱 III\ 🗉 🔕
	p Figure Tag Search Gallery Key-Value Search Admin	Search:	🔍 👤 Anna Ham
	Edit group Mame Demo_Group Description Demo_Group For demo and documentation purposes Fields marked in red are mandatory. Fields marked in red are mandatory. Permission: Private Read-Oniy Read-Annotate Read-Write Porners: Anna Hannacher (ansla 100) Bumers: Anna Hannacher	Cancel	

More information about the OMERO permission system and the group management can be found on the official support website: https://docs. openmicroscopy.org/omero/5.6.3/sysadmins/server-permissions.html and https://downloads.openmicroscopy.org/help/pdfs/group-owner.pdf.

Related articles

0

- OMERO upload (OMERO.insight)
 OMERO download
- OMERO search
- OMERO web linkOMERO data sharing