


Negative & Positive Stain (Waste Disposal)

Before work:

Where to find what and how to set up your Workspace...

1. "Waste Disposal-Kit" (under the fume hood)
2. staining solutions and stock solution at 4°C (the fridge right at the entrance)

Important: Only work within the labeled area on the appropriate surface in the designated workspace. If additional space is needed, kindly inform us, and we will prepare an alternative area accordingly

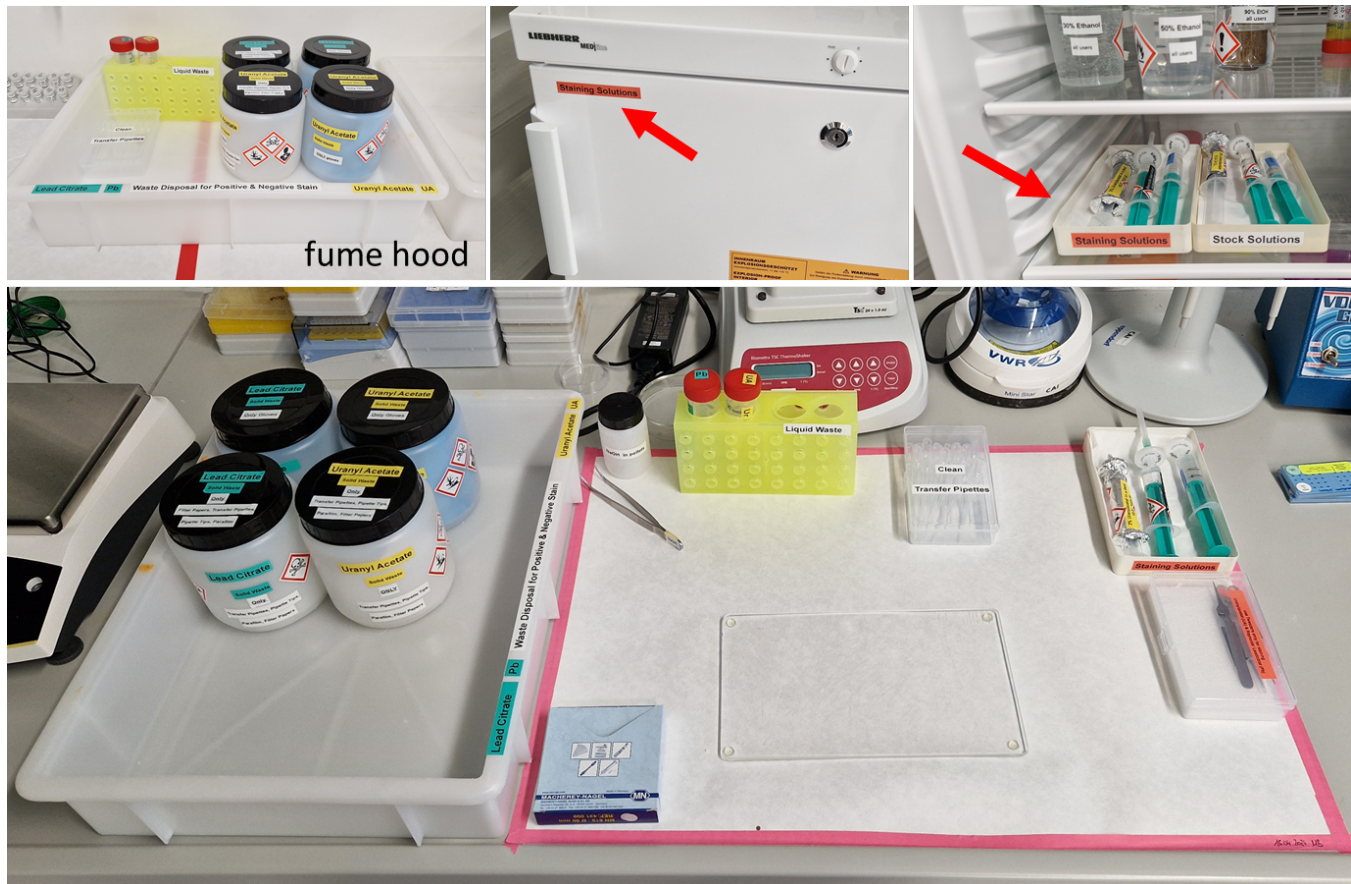


After work:

The procedure for cleaning up Uranyl Acetate is as follows:

(Lead citrate is cleaned up in the same manner, utilizing the corresponding containers in the waste disposal kit.)

1) Uranyl Acetate (UA) disposal

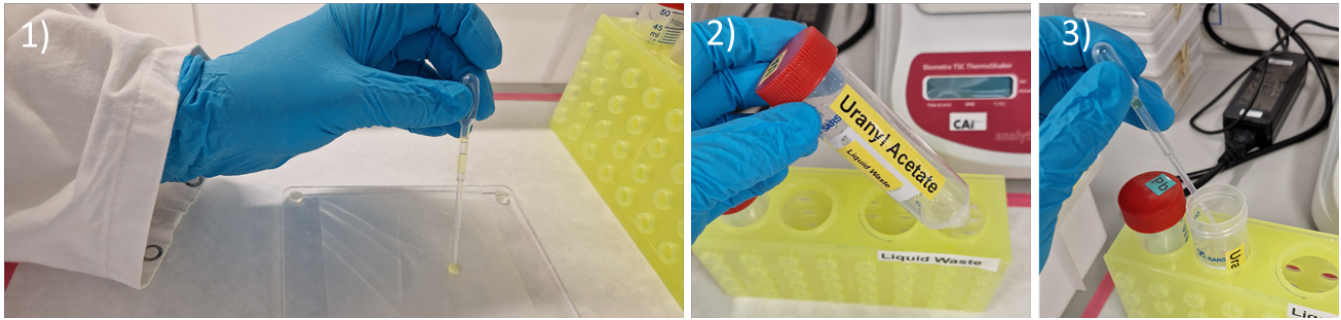

2) UA-Contaminated water droplets:

